

QUICK START GUIDE



Moticam T2

Moticam T2 Quick Start Guide

1. Remove the dust cap from the lens
2. Mount the Moticam T2 onto a C-mount adapter attached to your microscope
3. Plug in the power adapter into the Moticam T2
4. Press and hold the power button for a few seconds, wait until the Motic logo appears
5. Once it has finished loading, tap on the **MotiConnect T** app as shown below.



6. Now you should be able to see a live image on the screen of your Moticam T2




Connecting Your Device To The Moticom T2 Through Wi-Fi


1. Tap on the “Settings” button at the bottom



2. Make sure Wi-Fi is turned ON and select your network on the right, enter your Wi-Fi password (for protected network) then press “Connect”.

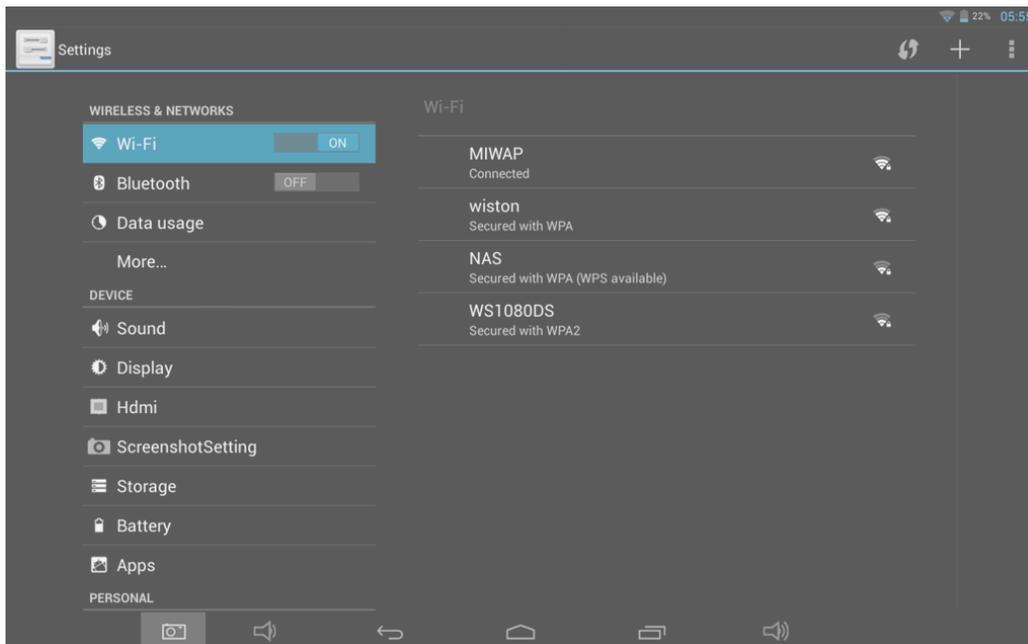

3. Go back to Moticonnect T app and press on the MoticHub button

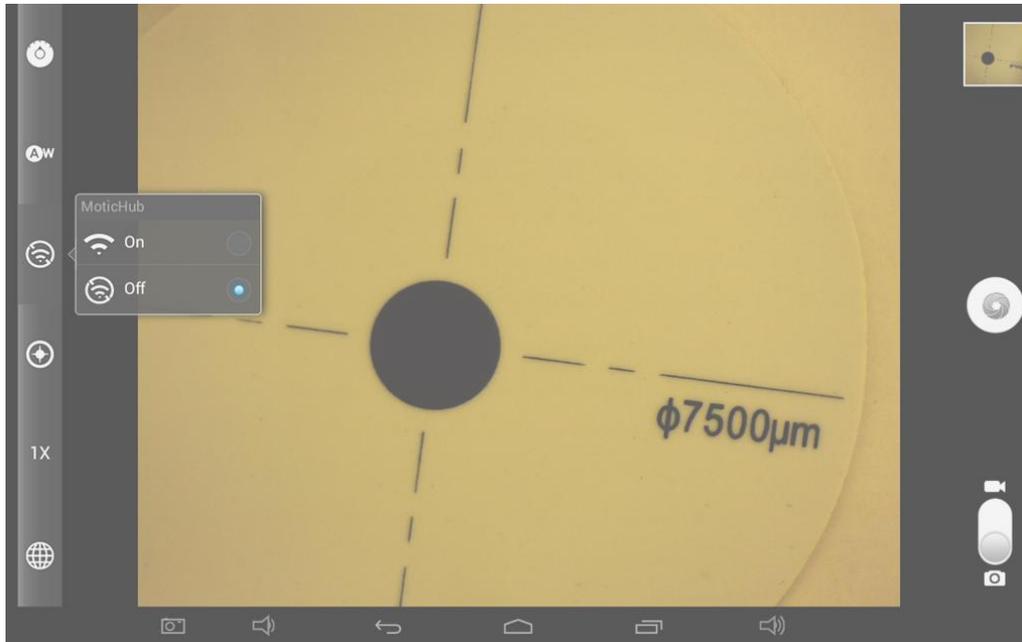

4. Select "On" to enable MoticHub on the Moticam T2. Use the IP Address in the dialog window displayed on the screen to connect any of your devices to the Moticam T2.



You can also check out the "Help" page for more details about the MotiConnect T app.

